

91600100 Rev D

# MERSCOPE<sup>®</sup> Vizualizer User Guide

*Desktop Version*

vizgen<sup>®</sup>

## NOTICES

Document number	91600100
Document revision	Rev D
Revision date	July 2024

### Legal notices

ALL PRODUCT(S) AND SERVICES DESCRIBED HEREIN ARE INTENDED FOR RESEARCH USE ONLY, AND NOT FOR USE IN DIAGNOSTIC OR TREATMENT PROCEDURES OR FOR ANY OTHER USE.

FAILURE TO COMPLETELY READ AND EXPLICITLY FOLLOW ALL OF THE INSTRUCTIONS CONTAINED HEREIN MAY RESULT IN DAMAGE TO THE PRODUCT(S), INJURY TO PERSONS, INCLUDING TO USERS OR OTHERS, AND DAMAGE TO OTHER PROPERTY. VIZGEN, INC. (“Vizgen”) DOES NOT ASSUME ANY LIABILITY ARISING OUT OF THE IMPROPER USE OF THE PRODUCT(S) DESCRIBED HEREIN (INCLUDING PARTS THEREOF OR SOFTWARE) OR ANY USE OF SUCH PRODUCT(S) OUTSIDE THE SCOPE OF THE EXPRESS WRITTEN LICENSES OR PERMISSIONS GRANTED BY VIZGEN IN CONNECTION WITH CUSTOMER’S ACQUISITION OF SUCH PRODUCT(S).

Nothing contained herein shall constitute any warranty, express or implied, as to the performance of any product(s) described herein. Any and all warranties applicable to any product(s) and software are set forth in the applicable terms and conditions of sale accompanying the purchase of such product(s). Vizgen provides no warranty and hereby disclaims any and all warranties as to the use of any third-party product(s) or protocols described herein. The use of product(s) described herein is subject to certain restrictions as set forth in the applicable terms and conditions of sale accompanying the purchase of such product(s).

Updates to product(s) and software may be required to enable customers to use product(s). If you fail to update the product(s) or software when provided by Vizgen, any warranty thereon shall immediately terminate. In the event of a product failure resulting from an update, such failed product will be subject to the warranty set forth in the terms and conditions agreed to in connection with the purchase of such product, if any, only if such product is covered by such warranty at the time of such failure. Vizgen is not required to repair, replace, or take any other remedial action with respect to any product not covered under such warranty.

Vizgen<sup>®</sup>, MERSCOPE<sup>®</sup> and MERSCOPE Ultra<sup>™</sup> are trademarks of Vizgen. All rights in the trademarks, service marks, trade dress, logos and copyrights are owned by Vizgen, Inc. and fully reserved. All other brands and names contained herein are the property of their respective owners.

This document and its contents are proprietary to Vizgen, Inc. and its affiliates, and are intended solely for the contractual use of its customer in connection with the use of the product(s) described herein and for no other purpose. This document and its contents shall not be used or distributed for any other purpose and/or otherwise communicated, disclosed, or reproduced in any way without the prior written consent of Vizgen. Vizgen does not convey any license under any patent, trademark, copyright, or common-law rights, nor similar rights of any third parties, by this document. The instructions in this document must be strictly and explicitly followed by qualified and properly trained personnel in order to ensure the proper and safe use of the product(s) described herein. All of the contents of this document must be fully read and understood prior to using such product(s).

© 2024 Vizgen, Inc. All rights reserved.

## SUPPORT

### Contact information

[support@vizgen.com](mailto:support@vizgen.com)

61 Moulton Street

West Cambridge Science Park Cambridge,

MA 02138

USA

Other Vizgen references are available online at <https://vizgen.com/>

## VERSION CHANGE NOTES

Summary of substantive changes from Rev C to Rev D

Brief description of change	Page(s)
Updated to accommodate compatibility with MERSCOPE Ultra Software (v234) and output file (.VZG2).	Throughout

## TABLE OF CONTENTS

<b>INTRODUCTION</b> .....	<b>7</b>
MERFISH Technology.....	7
The Vizgen MERSCOPE Platform Solution.....	7
The MERSCOPE Workflow.....	8
Broad Application.....	9
<b>VIZUALIZER SOFTWARE USER INTERFACE OVERVIEW</b> .....	<b>10</b>
<b>QUICK START ORIENTATION</b> .....	<b>11</b>
Installation.....	11
Create a First Experiment.....	11
Generate and Import Custom Cell Type Annotations.....	12
Open Custom Cell Clustering Analysis.....	13
<b>SYSTEM RECOMMENDATIONS</b> .....	<b>15</b>
<b>MENU BAR</b> .....	<b>16</b>
Software Versions.....	16
File > Assemble Experiment.....	16
File > Import Data.....	17
File > Import Data > Metadata.....	17
File > Import Data > Geometry > [Name of ROI Feature Set].....	19
File > Export Data.....	20
File > Export Data > Metadata.....	20
File > Export Data > Geometry > [Name of ROI Feature Set].....	21
Import/Export Leiden Colors From/To Scanpy.....	21
File > Open Experiment.....	21
File > Save Experiment.....	21
Edit > Manage Image Channels.....	22
Edit > Manage ROI Sets.....	22
Edit > Save Custom [Feature Set] Groups as New Copy.....	22
View.....	23
Help.....	24
Help > Archive Logs.....	24
<b>SPATIAL VIEW WINDOW</b> .....	<b>25</b>
Tooltips.....	25
Navigation and Views.....	25
Zoom and Scale.....	26
Information Panel.....	26
Export Image.....	27
Spatial Settings.....	27
Additional Settings (Under Spatial Settings Menu).....	28

<b>DETAILS AND SELECTED ITEMS PANELS.....</b>	<b>30</b>
Details Panel.....	30
Selected Items Panel - Cells.....	30
Selected Items Panel - ROI .....	31
<b>SPATIAL FEATURES PANEL – CELLS.....</b>	<b>32</b>
Select the Cells Feature Set.....	32
Cells .....	32
Active Category .....	33
Selection Toolbar (Cells).....	33
Cell Groups .....	35
Checked Cell Groups.....	37
Unchecked Cell Groups .....	37
Cell Group Settings .....	37
Cell Color Options.....	38
<b>SPATIAL FEATURES PANEL - REGIONS OF INTEREST.....</b>	<b>39</b>
Select an ROI Feature Set.....	39
Selection Toolbar (ROIs) .....	39
ROI Construction and ROI Groups .....	40
Checked ROI Groups.....	41
Unchecked ROI Groups .....	41
ROI Group Settings .....	41
ROI Color Options.....	42
<b>SCATTERPLOT .....</b>	<b>43</b>
Translate and Zoom.....	43
Select Array .....	43
Export Image .....	43
Selection Toolbar (Cells or ROI).....	44
<b>GENES.....</b>	<b>45</b>
Checked Genes .....	45
Unchecked Genes.....	45
Gene Color Settings .....	46
Gene Settings .....	46
Gene Highlighting.....	47
Gene Groups.....	48
Search for Genes .....	49
Gene Group Settings .....	49
<b>HEATMAP .....</b>	<b>50</b>
Navigation and Selection.....	50
Zoom and Scale .....	50
Export .....	50
Heatmap Toolbar.....	51
Heatmap Settings .....	52
Heatmap Highlighting.....	53

## INTRODUCTION

### MERFISH Technology

Multiplexed error-robust fluorescence in situ hybridization (MERFISH) is a spatially resolved single-cell transcriptome profiling technology. MERFISH combines the power of single-cell transcriptomics with spatial biology by directly visualizing and counting RNA transcripts from 100s to >10,000 genes in cells or tissue sections.

MERFISH advances the power of single-molecule fluorescence in situ hybridization (smFISH) – the gold standard of RNA quantification – with error robust barcoding, combinatorial labeling, and sequential imaging to greatly expand the multiplexing capacity. MERFISH enables researchers to map the molecular, cellular, and functional composition of biological systems with spatial context.

### The Vizgen MERSCOPE<sup>®</sup> and MERSCOPE Ultra<sup>™</sup> Platform Solutions

The Vizgen MERSCOPE Platform Solution is comprised of the Vizgen Gene Panel Design Portal, MERSCOPE reagent kits, the MERSCOPE Instrument, the MERSCOPE Analysis Computer, and the MERSCOPE Vizualizer software to streamline the acquisition of high quality MERFISH data and data interpretation.

The MERSCOPE Ultra Platform is Vizgen's latest generation spatial imaging platform. MERSCOPE Ultra offers the same high-quality, high-resolution MERFISH data as MERSCOPE but on a larger scale with an expansive imaging area up to 3.0cm<sup>2</sup>. MERSCOPE Ultra is comprised of the MERSCOPE Ultra Instrument and MERSCOPE Ultra Analysis PC and its workflow is also supported by the Vizgen Gene Panel Design Portal and MERSCOPE Vizualizer software.

Reagent kits facilitate sample preparation. The automated MERSCOPE and MERSCOPE Ultra Instruments integrate high-resolution imaging and fluidics to automatically acquire a full MERFISH dataset. The MERSCOPE Vizualizer software automates image processing and offers interactive visualization tools to explore MERFISH data.



### The MERSCOPE Workflow

The MERSCOPE Workflow begins with either designing a MERSCOPE Custom Gene Panel or ordering a MERSCOPE Predesigned Gene Panel.

**Design a custom gene panel.** The intuitive Vizgen Gene Panel Design Portal makes recommendations on your custom gene panel for optimal performance in a MERFISH measurement. For example, the portal highlights genes that might cause optical crowding or cannot be accurately targeted. Vizgen is currently offering custom gene panels for up to 140 genes, up to 300 genes, up to 500 genes, and up to 1000 genes.

**Order your MERSCOPE Gene Panel.** MERSCOPE Custom Gene Panels and MERSCOPE Predesigned Gene Panels are delivered in a ready-to-use format. Once the gene panel is finalized, users will be able to download their panel-specific MERSCOPE Codebook to the MERSCOPE Instrument.

**Prepare your sample.** The Vizgen sample preparation user guides take users through each step of the MERSCOPE sample preparation process.

**Load and run the MERSCOPE Instrument.** The MERSCOPE Slide is assembled into the MERSCOPE Flow Chamber and then loaded into the instrument along with a MERSCOPE Imaging Cartridge. Users define regions of interest on the MERSCOPE Slide and initiate the fully automated instrument run.

**Data Processing and Visualization.** The MERSCOPE Instrument Software (in combination with the MERSCOPE Analysis Computer) automatically processes the raw images to output spatial genomics measurements in a format ready for immediate downstream analysis. The output includes the list of all detected transcripts and their spatial locations in three dimensions (CSV files), mosaic images (TIFF), experiment metadata (JSON), output from the cell segmentation analysis: transcripts per cell matrix (CSV), cell metadata (CSV), cell boundaries (HDF5 [MERSCOPE Instrument Software v231 or

earlier] or PARQUET [v232 or later]), and a binary for use with the MERSCOPE Vizualizer software. The MERSCOPE Platform Solution includes the MERSCOPE Vizualizer software for visualizing and analyzing data. The output files are also compatible with open-source tools for single-cell and spatial analysis.

### Broad Application

The MERSCOPE Platform Solution has broad application in both fundamental biology and medicine – from basic science, to drug discovery, to clinical pathology. Find out more at

<https://vizgen.com/applications/>.



Oncology



Immunology



Neuroscience



Infectious Disease




Developmental Biology &  
Regenerative Medicine

## VIZUALIZER SOFTWARE USER INTERFACE OVERVIEW

1 Main menu bar	<p><b>File</b> Open experiment, recent, import data, export data, save experiment, assemble experiment</p> <p><b>Edit</b> Manage image channels and region of interest (ROI) sets, save a copy of custom feature set groups</p> <p><b>View</b> Control the visibility of the floating panels surrounding the spatial view window</p> <p><b>Help</b> Version information, about Qt (the C++ framework used for application development), open the report folder, archive logs</p>
2 Spatial view window	Tissue visualization, view the MERSCOPE Instrument output data
3 Heatmap	A two-dimensional view of cells by gene count data
4 Spatial features	Manage the active category, cell groups, and ROIs. Each feature set (cells, ROI sets) is addressed individually
5 Genes	View, select, and create gene groups
6 Scatterplot	Display two-dimensional arrays, such as cells in dimensionality reduced gene expression space (e.g., UMAP/tSNE*) or other user- imported two-dimensional arrays (e.g., cell volume versus transcript count)
7 Details	Information on a selected gene, cell, or ROI
8 Selected items	A list of selected cells and ROIs, and create cell and ROI groups

\*UMAP, uniform manifold approximation and projection; tSNE, t-distributed stochastic neighbor embedding.

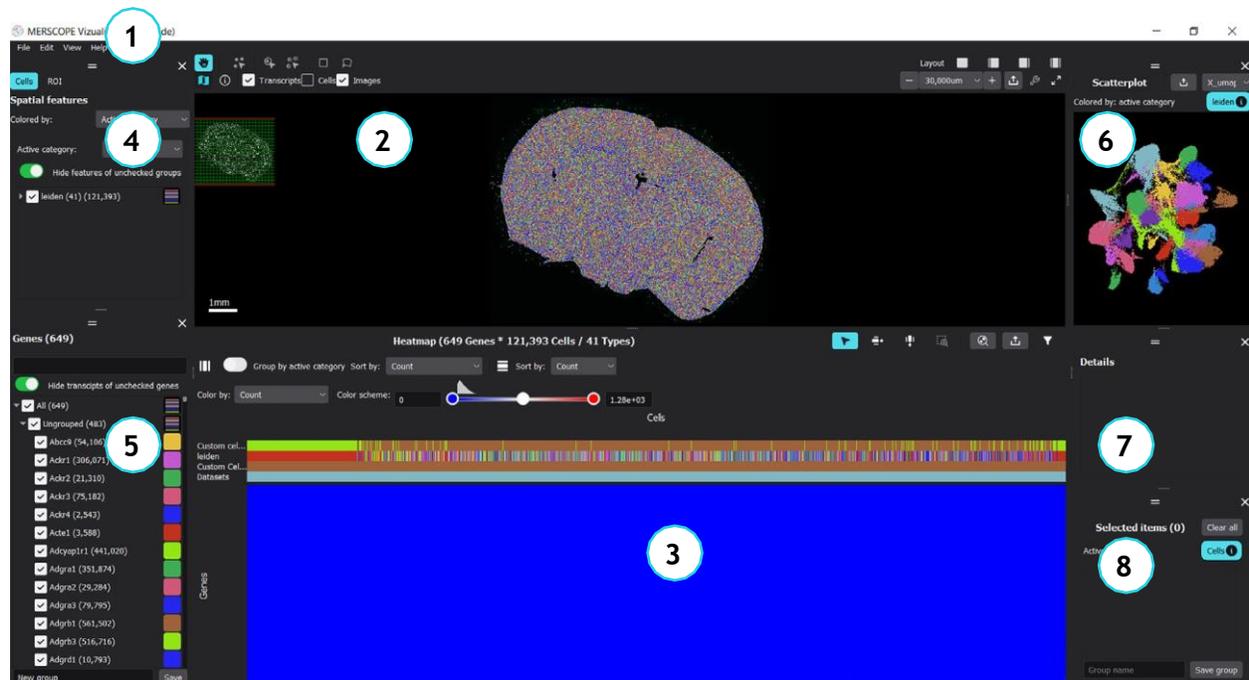

## QUICK START ORIENTATION

The MERSCOPE Vizualizer software enables interactive visualization and exploration of all components of a MERFISH measurement acquired on the MERSCOPE Instrument, including the detected transcripts and the segmented cell boundaries overlaid on the underlying images. The MERSCOPE Vizualizer uses data in the .VZG format, which is a compressed representation of the MERSCOPE Instrument output data optimized for efficient visualization.

Note: For instruments running control software version 234 and above, the .VZG format has been replaced with .VZG2. This new format includes more efficient file packing and stores fewer mosaic tiff images for a reduced, more manageable file size. Datasets from .VZG and .VZG2 can be combined in a single experiment within Vizualizer Software.

- *The quick start orientation narrative is intended to help users quickly understand basic functions and get started with the MERSCOPE Vizualizer software. Links to more detailed descriptions of software operations and additional features are included in the narrative.*

### Installation

- Install the MERSCOPE Vizualizer software using the installer executable and follow the installation prompts.
- Installation requires registering your installation with Vizgen. To register, follow the prompts, insert an email address, and enter the activation code that is shown on the registration landing page and emailed to you as a copy. Vizgen will use the provided email address to send notifications about new MERSCOPE Vizualizer software releases. Check [Software Versions](#) for any potential incompatibility between experiments assembled in different versions.
- *DO NOT install the MERSCOPE Vizualizer software on the MERSCOPE Instrument or the MERSCOPE Analysis Computer.*

### Create a First Experiment

- An experiment in the MERSCOPE Vizualizer software is a collection of datasets, where each dataset is a ROI defined on the MERSCOPE Instrument software during data acquisition. An experiment can contain one or more datasets that use the same MERSCOPE Gene Panel.
- A dataset .VZG/.VZG2 file can be found in the merfish\_output directory on the Z: drive of the MERSCOPE Instrument. Copy the .VZG/.VZG2 file to your computer either over the network or via a USB drive. → Open the MERSCOPE Vizualizer software and assemble the experiment by selecting File > Assemble experiment from the [MENU BAR](#).
- Enter a name for the new experiment in the Experiment name text field. Click the folder next to Select datasets to open a file browser and select the .VZG/.VZG2 file(s) copied from the MERSCOPE Instrument. Select an output location and click Build to assemble the experiment.

- The MERSCOPE Vizualizer software will unpack the .VZG/copy the .VZG2 file(s) into the specified experiment directory. Once the experiment is created, the user interface will populate with the experiment data.
- When creating an experiment from .VZG2 files, users are given the option to “Skip copying datasets” to save on disk space. When selected, this feature will create a hard link directly to the source .VZG2 file without creating a copy of the data. The hard link appears in the Experiment folder as a typical file but is only a pointer to the source data. As such, deleting the source file in its original location alone will not free up disk space. Deleting both the source file and hard link will recover disk space. When this feature is not selected, the software will duplicate the data from the source .VZG2 into the Experiments folder doubling the disk space requirement for the given experiment.
- Initially, the center z-plane of transcripts, cells, and images are displayed in the [SPATIAL VIEW WINDOW](#) (refer to [Spatial Settings](#) for more information on z-planes). All transcripts can be toggled on and off in the [Spatial Settings](#) menu (the tools icon above the top right of the spatial view window) and selecting All in the Z-plane menu for the transcripts layers. Cells and images can also be toggled on and off in this menu. All cell polygons can be visualized by selecting ‘All’ in Z-plane menu, while for images layer, only active z plane can be selected.  
  
NOTE: With .VZG2 files, transcripts from all z planes are represented in a single, stacked image and displayed as “All” always.
- To quickly visualize the measured expression of each gene, click on the name of a gene in the [GENES TREE](#) in the genes panel and use the up/down arrow keys to scroll through the list. When a gene is highlighted in the genes list, only the highlighted gene is displayed in color (the others are displayed gray) in the spatial view window (the transcripts layer must be checked). Press Esc to reset the selection.
- *Note that ROIs on the MERSCOPE Instrument are different from ROIs defined in the MERSCOPE Vizualizer software in the [SPATIAL FEATURES PANEL - REGIONS OF INTEREST](#).*

### Generate and Import Custom Cell Type Annotations

- The MERSCOPE Vizualizer software is able to visualize custom cell type annotations. To generate custom cell type annotations and populate the [SCATTERPLOT](#) with a UMAP representation of the data, the cell by gene matrix can be exported from the MERSCOPE Vizualizer software, loaded into custom analysis tools, exported with cell type annotations, and imported back into the MERSCOPE Vizualizer software with cell type annotations.  
MERSCOPE Vizualizer exports and imports cell annotations in the AnnData (HDF5), CSV, or PARQUET formats.
- To export an AnnData (.HDF5) file, select File > Export data > Metadata > [feature set] from the [MENU BAR](#). Select AnnData data format, navigate to the save location, and click Export.

## Open Custom Cell Clustering Analysis

- To execute this process, uploaded files must NOT have a `_clustered` suffix to the `.HDF5` file name (an error will occur during `Load AnnData`). If a file has previously been processed (clustered) and the file name changed to remove `_clustered`, errors may occur during processing.


→ A Vizgen template Jupyter Notebook is hosted on Google CoLab. To cluster the cells in the exported `AnnData` file in the Google CoLab notebook, navigate to <https://vizgen.com/merscope-clustering/> and click ACCESS THIS RESOURCE.

*Note: Google CoLab is a free service with a RAM limitation of approximately 13 GB, which may restrict the processing of larger datasets containing over 100,000 cells and 500 genes.*


→ Make a copy of the notebook in your account by selecting File > Save a copy in Drive.

→ Click on the folder icon (1) to the left of the browser window to open the Files tab.

→ When the Files tab is open, click on Upload to session storage (2) to upload a `.HDF5` file.

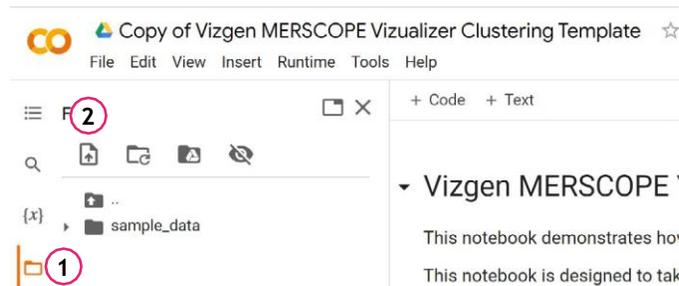


→ Click Refresh (3) to ensure the file is available and shown in the list.

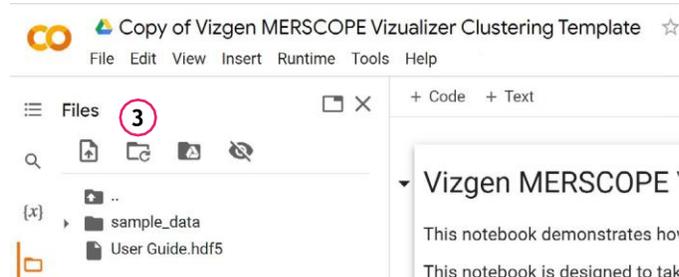

→ After the file has uploaded (note the progress wheel at the bottom of the Files tab), select Runtime > Run all (4) to execute all cells in the notebook.



- The notebook will run clustering analysis and save the clustered file with a `_clustered` suffix to the file name (it may take some time to appear). Right-click on the file name to download (5) the clustered AnnData file from the CoLab instance.

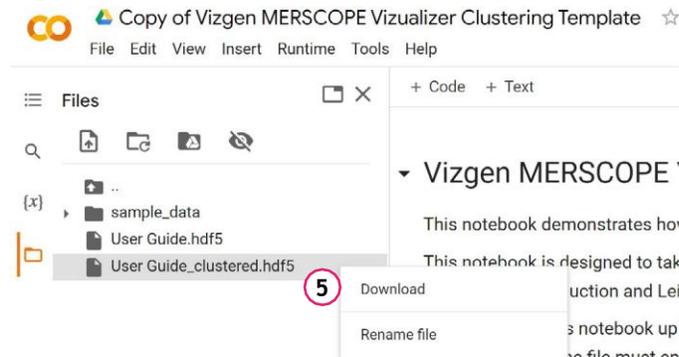

- Import the clustered AnnData file into the MERSCOPE Vizualizer by selecting File > Import data > Metadata > [feature set] from the [MENU BAR](#). Select AnnData data format, navigate to the AnnData file, and click Import.
- Both the clustering results and the two-dimensional representation of gene expression (e.g., UMAP) can be imported.
  - *There are many parameters that can be adjusted to obtain the most biologically relevant cell clustering parameters for a given dataset. For a new dataset, the cell filtering criteria and clustering algorithm parameters should be carefully considered.*
- After the clustering annotations and UMAP are imported, the [SCATTERPLOT](#) will display the imported UMAP and a new category corresponding to the new clustering annotations is available and selectable from the [SPATIAL FEATURES PANEL - CELLS](#). To color the cells using the clustering annotations, select the clustering annotations in the Active category menu, select Active category in the Color by menu, and ensure the cells layer is toggled on in the [Spatial Settings](#) menu.
- The gene expression for each cluster can be displayed in the [HEATMAP](#). Toggle on the Group by active category slider in the [Heatmap Toolbar](#).
- The UMAP can be colored by gene expression by selecting Count or Z-score (among others) from the Color by menu in the [SPATIAL FEATURES PANEL - CELLS](#).

## SYSTEM RECOMMENDATIONS

For optimal performance, the MERSCOPE Vizualizer software should be run on a system with the following minimum criteria:

- Processor: Intel Core i9
- RAM: 32 GB
- Graphics: NVIDIA GeForce RTX 3080-10 Gb
- SSD Capacity: 1 TB

The MERSCOPE Vizualizer software can be run on a system with lower capability, but may result in suboptimal performance, especially for large datasets.

macOS 12 (Monterey) is recommended, where applicable.

In the absence of an optimal graphics card, the MERSCOPE Vizualizer software can be run on a GPU-enabled AWS EC2 instance. Install NVIDIA drivers following the instructions under Option 4: NVIDIA gaming drivers (G5 and G4dn instances) at the following link:

- <https://docs.aws.amazon.com/AWSEC2/latest/WindowsGuide/install-nvidia-driver.html#nvidia-installation-options>

## MENU BAR

### Software Versions

After an experiment is assembled in version 2.4, it cannot be opened in earlier versions because the folder structure of the assembled experiments is different. Experiments created in version 2.3 are compatible with version 2.4. The MERSCOPE Vizualizer software version is found under the [Help](#) menu.

### File > Assemble Experiment

To assemble an experiment:

- File > Assemble experiment.
- Enter Experiment name (this is then shown in the top bar of the MERSCOPE Vizualizer software window).
- Click the Select output folder icon to save MERSCOPE Vizualizer metadata.
- Click the Select datasets folder icon to select datasets (.VZG/.VZG2 files).
- Click Build.
- *Multiple datasets can be assembled, but they must be from samples measured with the same MERSCOPE Gene Panel. To assemble multiple datasets, repeatedly click Select datasets and select the datasets for inclusion in the experiment. If datasets are in the same folder, use CTRL-SHIFT-select to simultaneously select multiple datasets.*
- *Assembling an experiment may take some time. A progress bar across the top of the spatial view window shows progress.*
- *The MERSCOPE Vizualizer software creates a new folder with the experiment name in the output folder.*
- *Only the transcripts and images layers are visible by default in the spatial view window when an experiment is loaded. Refer to [Spatial Settings](#) for more information.*

## File > Import Data

The MERSCOPE Vizualizer software supports the import of data to feature sets via:

- The open HDF5 format: <https://www.hdfgroup.org/solutions/hdf5/> (annotated data, AnnData).
- .CSV files or .PARQUET files (categories files and numeric categories files). A numeric categories file imports data to the scatterplot. Data in .CSV and .PARQUET files must be organized in the same way as exports from the MERSCOPE Vizualizer software.
- .CSV files of ROI geometry data (of ROI feature sets only - this option is not available for the cells feature set).

## File > Import Data > Metadata

After processing and analyzing annotated data, it can be imported into the MERSCOPE Vizualizer software with additional information (e.g., clusterization).

Each experiment contains two default cell categories:

- Datasets is an immutable cell category. This category contains all the cells in each dataset.
- Custom Cell Groups. When an experiment is first assembled, a single cell group exists in this category (unclustered), which contains all the cells in a dataset. Custom cell groups can be created within this category.

Each experiment contains one default ROI category:

- Custom ROI Groups. When an experiment is first assembled, a single ROI group exists. Custom ROI groups can be created within this category.

In addition to default cell/ROI categories, additional cell/ROI categories can be imported with annotated data (e.g., Leiden). Cell/ROI categories can be linked (nested) during import.

In an experiment assembled from multiple datasets (.VZG/.VZG2 files), an exported metadata file contains all cells/ROIs from all datasets in the experiments, combined. In an imported metadata file, the clustering applies to all cells/ROIs in the experiment.

### To import annotated data (default):

– *Advanced import options are on the following page. If annotated data has previously been imported, the advanced Categories menu automatically appears.*

→ File > Import data > Metadata > select the spatial feature set (cells or an ROI) to import data into.

- Select the data format (AnnData [.HDF5], .CSV, or .PARQUET), navigate to the file location, and click Import. A Categories dialog box appears.
- Check the categories for import.
- Data available for the scatterplot (e.g., UMAP, tSNE) are imported by default, and can be renamed in the dialog box (under two-dimensional arrays for scatterplot). If not required, uncheck the appropriate option. Multiple selections can be checked/unchecked and a menu in the scatterplot window controls the data being displayed.
- Click Import.
- *The imported cell categories appear in the active category selection menu in the cells feature set in the spatial features panel and the scatterplot is populated.*
- *If a spatial feature set is empty (no features) an error message appears.*

**To import annotated data (advanced):**

- File > Import data > Metadata > select the spatial feature set (cells or ROI) to import data into.
- Select the data format of the file (AnnData [.HDF5], .CSV, or .PARQUET), navigate to the file location, and click Import. A Categories dialog box appears.
- Click the Show advanced settings button (if available - if annotated data has previously been imported, the advanced Categories menu automatically appears).
- Check the categories for import.
- Check a cell/ROI category in the Existing experiment metadata list to import annotated data for that category.
- *The default cell/ROI categories (datasets and custom cell/ROI groups) are already selected in the existing experiment metadata list.*
- Cell/ROI categories can be linked. Select a subcategory to nest it under a category.
- *Linking imported cell/ROI categories may be useful when there is a connection between them - e.g., coarse and fine clusters.*
- Cell/ROI categories can be renamed. Click on the name and enter the new name. Each category should have a unique name.
- *If a previously used name is entered, the MERSCOPE Vizualizer software automatically renames the category to be imported to give it a unique name.*

- *Default cell categories (datasets and custom cell groups) cannot be renamed or linked. Default ROI categories can be removed only if the entire ROI feature set is removed.*
- Data available for the scatterplot (e.g., UMAP, tSNE) are imported by default, and can be renamed in the dialog box (under two-dimensional arrays for scatterplot). If not required, uncheck the appropriate option. Multiple selections can be checked/unchecked and a menu in the scatterplot window controls the data being displayed.
- Click Import.
- *The imported cell/ROI categories appear in the active category menu in the feature set's spatial features panel and the scatterplot is populated (as applicable).*
- *If a ROI feature set is empty (no features) an error message appears.*

**File > Import Data > Geometry > [Name of ROI Feature Set]**

To import the coordinates of ROI polygon vertices (coordinates should be in well-known text [WKT] format):

- File > Import data > Geometry > [Name of ROI feature set].
- Navigate to the .CSV files for import.
- Click Open.
- *The ROI groups will appear in the spatial features panel and the spatial view window.*

## File > Export Data

The MERSCOPE Vizualizer software supports the export of data from feature sets to:

- The open HDF5 format: <https://www.hdfgroup.org/solutions/hdf5/> (annotated data, AnnData).
- .CSV files or .PARQUET files (expression matrix files, categories files, and numeric categories files for cells and ROI feature sets; and feature metadata files for ROI feature sets). The feature metadata files contain feature volume/area and coordinates.
- .CSV files of ROI geometries (to ROI feature sets only - this option is not available for the cells feature set).

## File > Export Data > Metadata

Exported annotated data can be processed and analyzed in third-party tools for analyzing single-cell gene expression data (e.g., Scanpy <https://scanpy.readthedocs.io/en/latest/#>).

In an experiment assembled from multiple datasets (.VZG/.VZG2 files), an exported metadata file contains all cells from all datasets in the experiments, combined.

### To export annotated data:

- File > Export data > Metadata > Select the spatial feature set (cells or ROI) to export data from.
- Select the format of the file (AnnData [.HDF5], CSV, or PARQUET).
- For export to .CSV and .PARQUET files: Check the files for export (expression matrix file, categories file, numeric categories file, and/or feature metadata file) and navigate to the location to save the files. Select the data type for the expression matrix file (count, z-score, normalized count, normalized z-score) from the menu.
- Confirm all save locations and file names are per preferences.
- Click Export.

### File > Export Data > Geometry > [Name of ROI Feature Set]

#### To export the coordinates of ROI polygon vertices:

- File > Export data > Geometry > [Name of ROI feature set].
- Navigate to the location to save the .CSV files.
- Click Save.

### Import/Export Leiden Colors From/To Scanpy

If a color scheme is applied in Scanpy (e.g., Leiden), the color scheme will be imported with annotated data file and displayed in the MERSCOPE Vizualizer software. Any changes made to the colors in the MERSCOPE Vizualizer software will be exported in the AnnData file and may be imported into Scanpy.

### File > Open Experiment

To open an experiment:

- File > Open experiment.
- Select the folder with the experiment data.

It is also possible to open a recent experiment via Recent > [name of experiment].

- *The MERSCOPE Vizualizer software opens the data folder and displays the experiment. If the experiment has annotated data, the software also loads all annotations.*

### File > Save Experiment

Assembled experiments are auto-saved after any changes are made in the application. All imported data and data created in the MERSCOPE Vizualizer software (custom cell groups, ROIs, gene groups, custom colors of cell/ROI/gene groups) are saved.

To manually save an experiment at any time:

- File > Save experiment.
- *All imported data and data created in the MERSCOPE Vizualizer software are automatically saved on closing the application.*

### Edit > Manage Image Channels

Manage the image channels available in the experiment by unchecking/checking the boxes next to each channel. DAPI and PolyT channels are included in all experiments. Other imaging channels (up to 16) are available for import based on the use of auxiliary cell boundary stains/other auxiliary stains during data acquisition. Selected image channels are listed in the [Spatial Settings](#) menu.

### Edit > Manage ROI Sets

ROIs can be defined. There is one empty ROI feature set by default (named ROI). ROIs can be grouped in each ROI feature set.

#### To manage ROI feature sets:

- Edit > Manage ROI sets.
- Click Add feature set to add a new ROI set.
  - *A new ROI set with a default name appears in the dialog box. The default name can be changed in the dialog box. The name of the default ROI set (ROI) can also be changed from the dialog box. The ROI set name should be unique.*
- Rename the ROI set or accept the default name and click OK.
  - *Multiple ROI sets can be added from the same dialog box.*
  - *A new ROI set appears in the spatial features panel.*
  - *ROI sets can be deleted or renamed from the manage ROI sets dialog box.*

### Edit > Save Custom [Feature Set] Groups as New Copy

#### To save custom cells/ROI groups as a new copy:

- Edit > Save Custom Cells groups as new copy
- Insert a name for the copy and click OK.
  - *The copied and renamed custom cells/ROI groups can be selected in the spatial features panel from the [Active Category](#) menu. The copied group is immutable (only group colors can be changed by the user).*


## View

Floating panels surround the spatial view window (refer to [VIZUALIZER SOFTWARE USER INTERFACE OVERVIEW](#) for the default layout).

- Scatterplot
  - Spatial features
  - FPS measuring (FPS = frames per second)
  - Genes
  - Details
  - Selected items
- All panels except FPS measuring are shown by default after launching the MERSCOPE Vizualizer software.

### To manage floating panels:

→ View > toggle the applicable panel on and off.

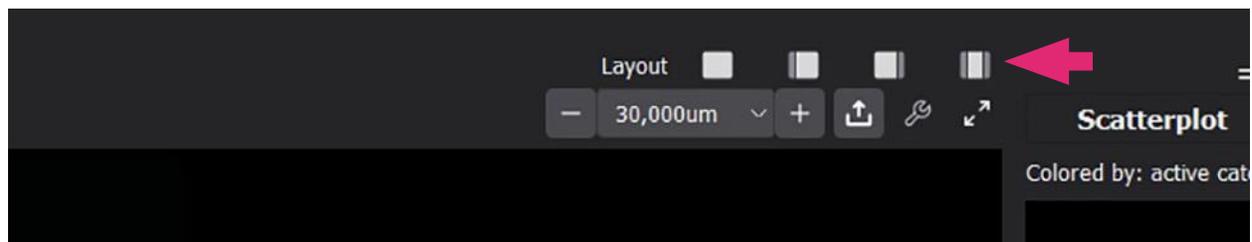

- To change the relative layout of the spatial view window and panels, toggle among the layout icons above the top right of the spatial view window.
- To move a panel, move the cursor toward the top of the panel until the hand cursor appears. Click and drag to move the panel. Blue highlighting appears when hovering over a new location.
- To create layered floating panels, move the cursor toward the top of the panel until the hand cursor appears. Pick up the panel, drag and drop it over an existing floating panel (blue highlighting appears when hovering over the new location). The layered panels can be selected from tabs at the bottom of the window. To return to a classic layout, toggle among the layout icons above the top right of the spatial view window.
- To hide a panel, click X in the top right corner or toggle off from the View menu.
- To resize a panel, hover over the bottom of a panel until the resize cursor appears. Click and drag to resize.

- When the FPS measuring panel is open, click Measure FPS and a number appears after some time (>20 is reasonable, >60 is optimal).
- *If FPS is <5, the user experience may be improved by upgrading the graphics capabilities of the computer.*

## Help

- Contact Vizgen help via [support@vizgen.com](mailto:support@vizgen.com).
- About > version information.
- About Qt (the C++ framework used for application development).
- Open the report folder.

## Help > Archive Logs

To create a .ZIP folder to share with [support@vizgen.com](mailto:support@vizgen.com):

- Help > Archive logs.
- Describe the problem in the text box.
- Attach a screenshot, if applicable.
- Click Save.

The .ZIP folder will contain (as applicable):

- The application version.
- Information from the manifests of the last loaded experiments.
- Information from the manifests of the last assembled datasets.
- Debug information about application function.
- Crash dump files.

## SPATIAL VIEW WINDOW

The MERSCOPE Vizualizer software has the capacity for two spatial view windows. The properties of each window are controlled independently. Each spatial view window displays a dataset in the experiment. If only one dataset is included during experiment assembly, only one spatial view window is visible by default.


To open the second spatial view window, drag the right border of the spatial view window to the left. In the [Spatial Settings](#) menu of the new empty spatial view window, select the current dataset from the dataset menu (which is None by default). When two spatial views windows are open, it is possible to visualize the same dataset in the same or different locations.

In the spatial view windows, users can work with existing experimental data (cells, transcripts, images) and create custom cell groups and ROIs. Cell selection and cell group formation or ROI selection and ROI group formation in the spatial view window and scatterplot are synchronized.

### Tooltips

→ Hover over a transcript, cell, or ROI to reveal a tooltip with information about the item.

### Navigation and Views

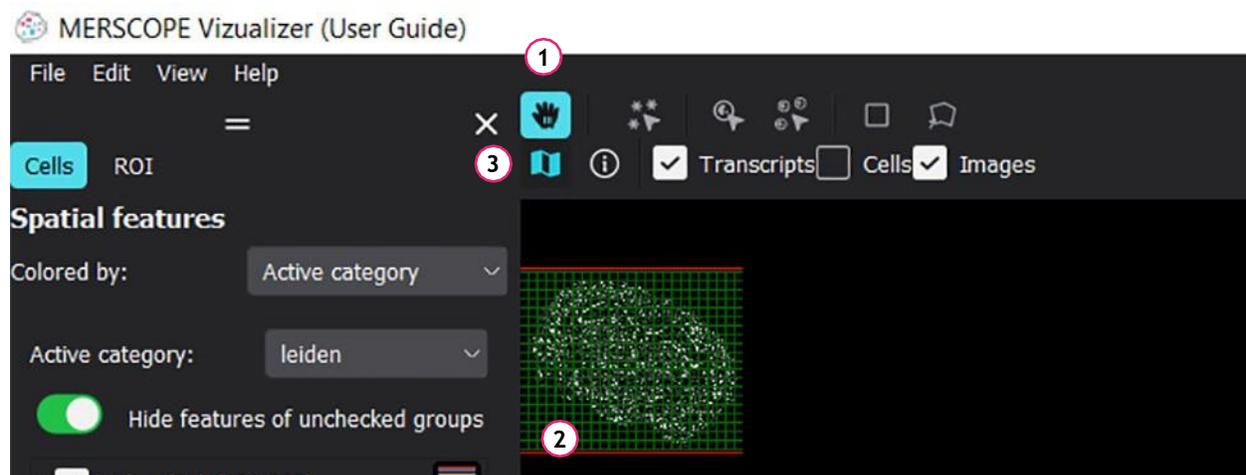


→ In the zoom/pan cursor mode (hand icon) (1), click and drag the hand cursor in the spatial view window and to move the content within the window.

→ There is a mini map (2) - a small scale preview of the spatial view window - in the top left corner of the window. Click and drag to move the mini map boundary (red box) to navigate the content of the spatial view window. This feature is particularly useful when zoomed in.


NOTE: This mini map feature is based on Images and changing the Image Channel Selection and brightness (under Spatial Settings menu) also changes the mini map. In Vizualizer versions 2.3 and below, the minimap is based on density.

→ Toggle the mini map on and off via the map icon above the top left of the spatial view window (3).

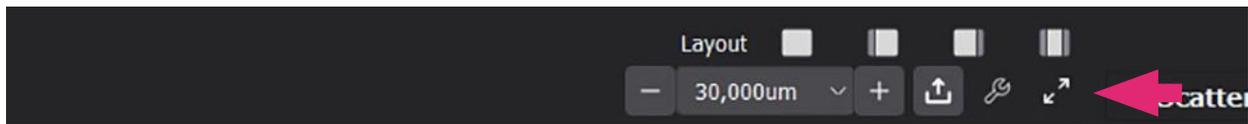

→ Change the content of the spatial view window to full screen and back by toggling the expand icon above the top right of the window.


## Zoom and Scale

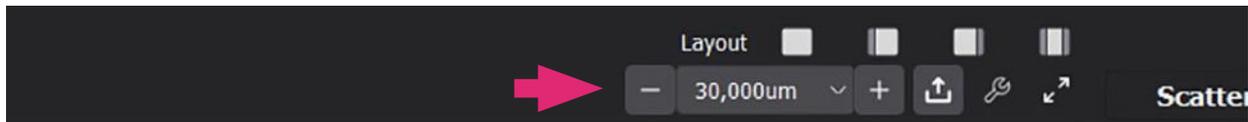

**To zoom in/out of the content in the spatial view window:**


- Rotate the mouse wheel.
- Click the +/- buttons to either side of the scale box above the top right of the window.
- Type in an exact zoom level in microns in the scale box. The value corresponds to the width of the spatial view window. A scale bar is always present in the bottom left of the window.
- Select a zoom level in microns from the menu in the scale box.

## Information Panel

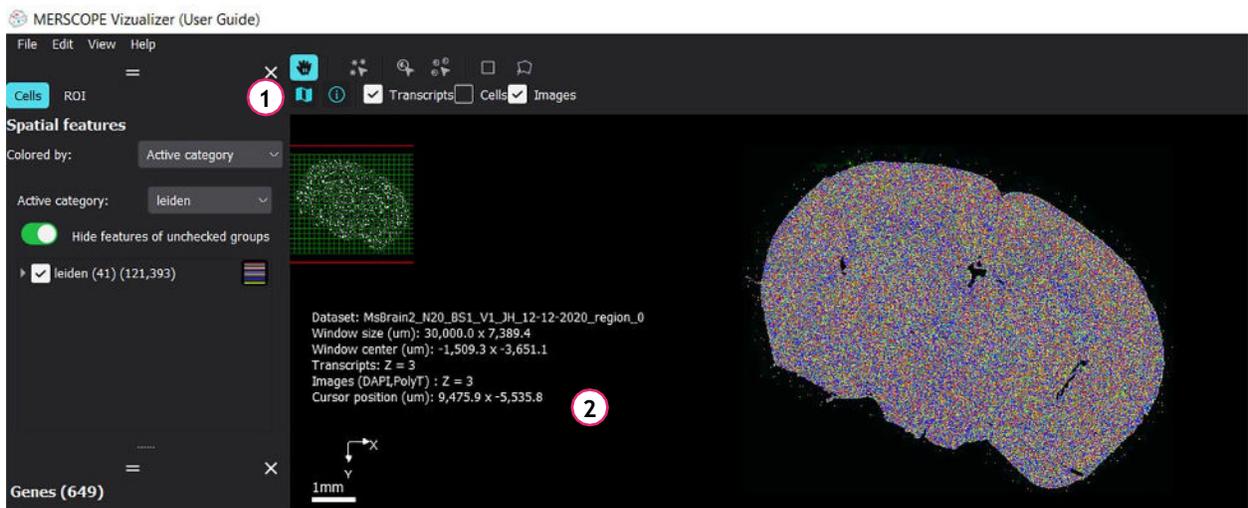


- Toggle a summary of properties of the data displayed in the window on and off via the information icon above the top left of the spatial view window (1).
- The information panel is displayed in the bottom left of the spatial view window (2), and includes the coordinates of the cursor in the spatial view window.

## Export Image

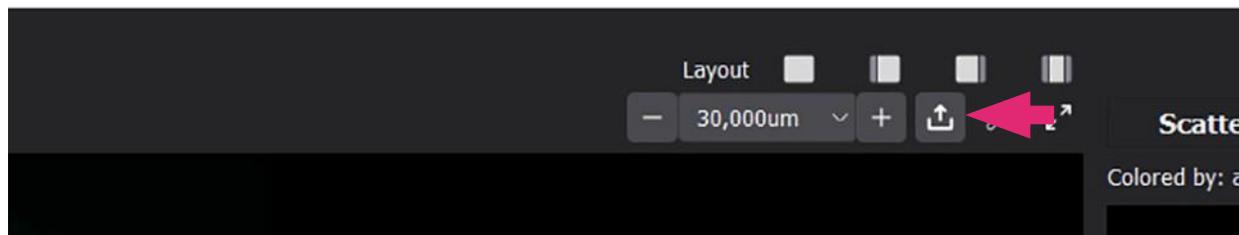

- Click the export icon above the top right of the spatial view window to export a .png file of the spatial view window. A .txt file with the same file name with details of the data in the image file is also created.
- Changing the resolution factor of the .png file changes the resolution and size of the image file. The range of resolution factors depends on the rendering capabilities of the computer and the actual size of the exported spatial view window (a smaller window can be exported with a higher resolution factor).

## Spatial Settings

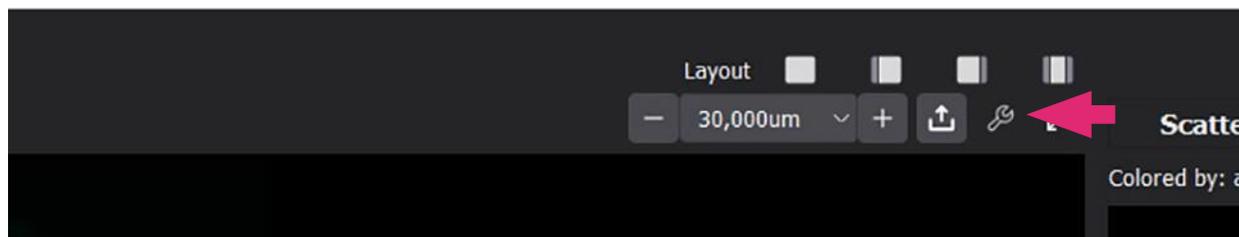

**To manage the properties of a spatial view window, toggle the spatial settings menu on and off via the tools icon above the top right of the spatial view window:**

- Select the dataset visible in each spatial view window in the Dataset menu.
- To choose the z-plane visible in the spatial view window, select from the Z-plane menu. The active (selected) z-plane or all z-planes can be selected for each layer. For .VZG2 files, only the active z-plane is available for Images layer.
- *Z-plane 0 includes the data at 1.5 microns from the MERSCOPE Slide. Z-planes 1, 2, 3, etc are at an additional fixed 1.5 microns away from the MERSCOPE Slide. The number of z-planes depends on the thickness of the sample defined during data acquisition.*

*NOTE: Transcripts are no longer divided by z planes for .VZG2 file format. Transcripts from all layers are represented in a single image.*

- Only the transcripts and images layers are visible in the spatial view window when an experiment is loaded. Additional layers are ROI feature sets, cells, and density. The density layer, however, is only available for VZG files.
- To toggle layers on and off in the spatial view window, check/uncheck the layer in the spatial settings menu or the shortlist of layers (transcripts, cells or the selected ROI feature set, images) above the top left of

→ **To manage the order of layers, click the Reorder layers button and a dialog box appears:**

- The transcripts layer is always on top (first on the list) so that the user can always select transcripts in the spatial view window.
- The active feature set is always listed second. Switching the active feature set in the spatial features panel automatically changes the layer list order. Selection tools always work on the active feature set.
- For the remaining layers, drag and drop a layer into a different location to reorder layers.

→ Each layer represents a particular type of spatial data. Cells and transcripts layers can be viewed across all z-planes or the active z-plane selected at the top of the spatial settings menu:

- **Transcripts.** Shows the checked gene transcripts as dots. A slider controls the intensity. A separate slider controls the size of the transcripts.
- **Cells.** Shows the polygons representing cells, identified during data processing. The slider controls the transparency.

NOTE: On distant zoom levels, Cell Selection by polygon/rectangle selects cells whether the current z-plane contains cell polygon or not. This is because cell selection is done using cell center coordinates. While on the close zoom levels, cell selection by polygon or rectangle selects cells whether current z-plane contains cell polygon or not, as in this case the selection is done using cell polygons.

- **ROI feature sets.** Shows the ROIs in the set. Each ROI feature set has its own layer. The slider controls the transparency. ROIs span all z-planes and an ROI has the same boundary for all z-planes.
- **Density.** Shows the density of checked gene transcripts. The slider controls the transparency. Density is available only for .VZG datasets.
- **Images.** Images of the image channels, if checked (e.g., DAPI and PolyT). Up to 16 image channels can be displayed at the same time ([Edit > Manage Image Channels](#)). The Images slider is used to set the overall transparency for the images. Individual channel sliders are used to set minimum and maximum contrast thresholds for each channel. Only one z-plane (active) can be selected in the image layer.

NOTE: All image channels will have the same number of z-planes for VZG2. Furthermore, images can have one z-plane (2D mode) and no z-plane switching in this case for VZG2.

### Additional Settings (Under Spatial Settings Menu)

- To define the coordinate at the center of the spatial view window, enter X and Y values.
- To rotate the content in the spatial view window, adjust the Rotate image setting. To rotate clockwise, enter a positive number. To rotate counterclockwise, enter a negative number.

- Click the Background color palette to change the background color of the window.
- *If image channels are selected and/or not transparent, the background color change is not visible until the images are deselected in the spatial settings menu or from the layers check boxes above the top left of the spatial view window.*

## DETAILS AND SELECTED ITEMS PANELS

### Details Panel

The details panel displays details of a selected gene, cell, or ROI, including a list of genes expressed in a selected cell/ROI. Data are organized under spatial and heatmap headings. The information can be copied to clipboard using the copy icon at the top right of the details panel.

### Selected Items Panel - Cells


- The selected items panel indicates the active feature set, i.e., the spatial feature set selected in the spatial features panel.
- When the cells feature set is active in the spatial features panel, all cells selected in the spatial features panel, the spatial view window, scatterplot, and/or heatmap are listed in the selected items panel.
- Cell indexes and identification numbers of the selected cells are also listed. The number of selected cells is listed in parentheses next to the panel title Selected items.
- Click Clear all on the selected items panel to clear the selection in all panels. The selected item list is empty if no cells are selected (when in the cells feature set).
- Switching between spatial feature sets (cells or ROI) clears the selected items list.
- Individual cells in the list can be selected by clicking the cell in the list and the cell's location is identified by crosshairs in the spatial view window, scatterplot, and heatmap. The location is identified in the heatmap as long as the Group by active category slider is toggled off in the [Heatmap Toolbar](#).
- Alternatively, copy the cell name (e.g., 300948112130576190679118935843036609088) to clipboard. In the selected items panel, right-click (or Ctrl+V) to paste the cell name, and then select the cell in the single-item list to identify (crosshairs) the cell in the spatial view window, scatterplot, and heatmap.
- Cell groups ([Cell Groups](#)) can be created from this panel.

### Selected Items Panel - ROI

- The selected items panel indicates the active feature set, i.e., the spatial feature set selected in the spatial features panel.
- When an ROI feature set is active in the spatial features panel, all ROIs selected in the spatial features panel, spatial view window, scatterplot, and/or heatmap are listed in the selected items panel.
- The number of selected ROIs in the group is listed in parentheses next to the panel title Selected items. Right-click on an ROI name to delete it.
- Click Clear all on the selected items panel to clear the selection in all panels. The selected item list is empty if no ROIs are selected (when in an ROI feature set).
- Switching between spatial feature sets (cells or ROI) clears the selected items list.
- Individual ROIs can be selected by clicking the ROI in the list and the center of the ROI is identified by crosshairs in the spatial view window, scatterplot, and heatmap. The location is identified in the heatmap as long as the Group by active category slider is toggled off in the [Heatmap Toolbar](#).
- ROI groups ([ROI Construction and ROI Groups](#)) can be created from this panel.

## SPATIAL FEATURES PANEL - CELLS

### Select the Cells Feature Set

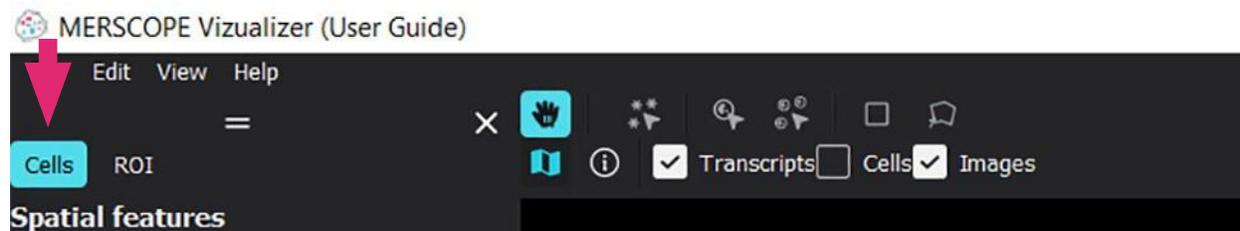

The cells feature set and ROI feature sets ([Select an ROI Feature Set](#)) are individually addressable by selection of a feature set at the top of the spatial features panel. The selected items panel also indicates the active feature set, i.e., the spatial feature set selected in the spatial features panel. The active feature set is always displayed in front of other feature sets in the spatial view window and selection tools are only applicable to the active feature set.

### Cells

The cell groups available in each cell category are listed in the cells feature set of the spatial features panel.

Each experiment contains two default cell categories:


- Datasets is an immutable cell category. This category contains all the cells in each dataset.
- Custom Cell Groups. When an experiment is first assembled, a single cell group exists in this category (unclustered), which contains all the cells in an experiment. Custom cell groups can be created within this category.

In addition to default cell categories, additional cell categories can be imported with annotated data (e.g., Leiden).

It is possible to make a copy of custom cell groups ([Edit > Save Custom \[Feature Set\] Groups as New Copy](#)).

## Active Category


Select an active category from the menu near the top of the spatial features panel. The menu contains the default categories and any imported cell categories. The data of the active category is displayed in the spatial view window, scatterplot, and heatmap.

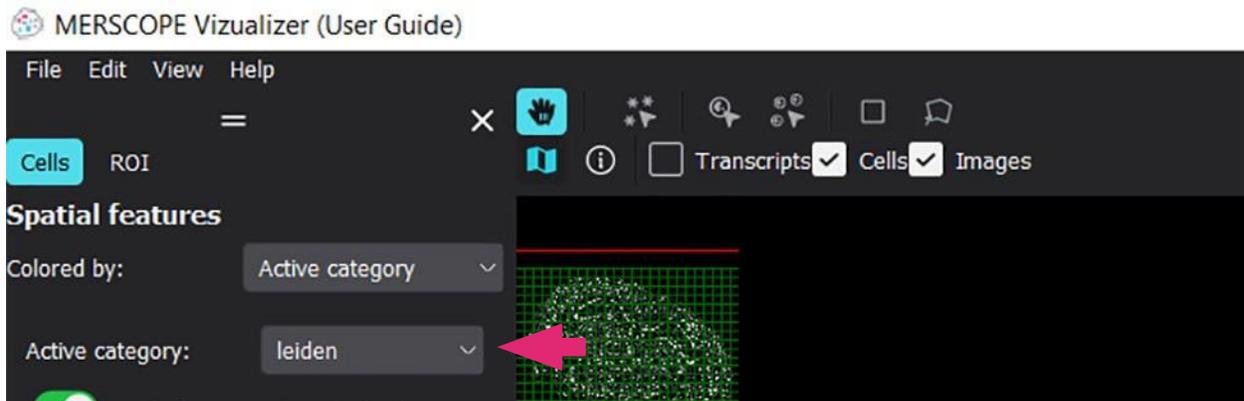

## Selection Toolbar (Cells)

The cells layer in the spatial view window must be turned on to select cells and define cell groups. To toggle the cells layer on and off, check/uncheck the cells layer in the [Spatial Settings](#) menu or in the shortlist of layers above the top left of the spatial view window (available when the cells feature set is selected).

A selection toolbar is available above the top left of the spatial view window; the selection tools also work in the scatterplot.

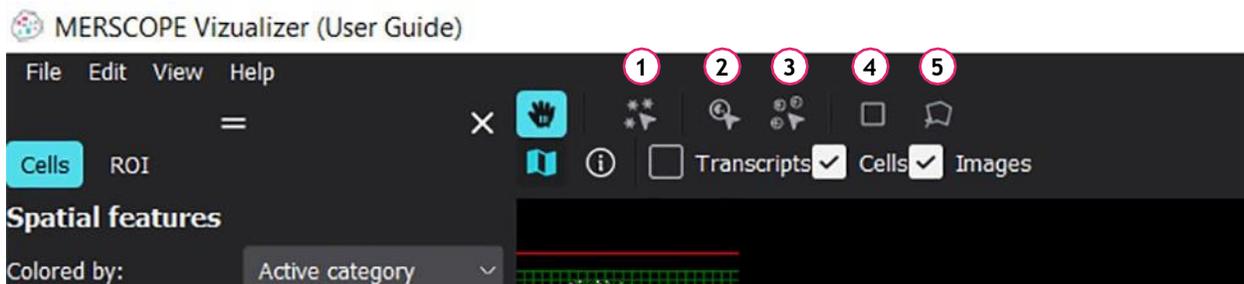

→ Multi-transcript selection (1) enables users to select all instances of a gene transcript at the same time. Click away from a selected transcript or press Esc and all [Checked Genes](#) transcripts reappear.

NOTE: For experiments created from .VZG2 files, multi-transcript selection works only with the closest Level of Details (LOD0).

→ Single-feature selection (Cells) (2) enables users to select a single cell. Click/click again to select/deselect an individual cell; click Clear all on the selected items panel, or press Esc to clear the selection. Select multiple cells by Ctrl-clicking single cells one-by-one. To deselect a cell from the selection, Ctrl-click in the spatial view window, scatterplot, or heatmap (as long as the Group by active category slider is toggled off in the [Heatmap Toolbar](#)).


- Multi-feature selection (Cells) (3) enables users to select all cells in a cell group. Click/click again to select/deselect a cell group; click Clear all on the selected items panel to clear the selection. Select multiple cell groups by Ctrl-clicking cell groups one-by-one.
- Rectangle selection (4). Drag and click to define a rectangle over the display to select cells at least partially within the rectangle. Press Ctrl and draw another rectangle to add cells to the previous selection. Draw another rectangle without Ctrl away from the original to reset the selection. Draw another rectangle within the first to select the subset of cells. Switch to the single-feature selection tool and click in a blank space in the spatial view window, click Clear all on the selected items panel, or press Esc to clear the selection.
- Polygonal lasso selection (5). Click each point to draw a polygon over the display. Close the polygon to select the cells at least partially within the polygon. Press Ctrl and draw another polygon to add cells to the previous selection. Draw another polygon without Ctrl away from the original to reset the selection. Draw another polygon within the first to select the subset of cells. Switch to the single-feature selection tool and click in a blank space in the spatial view window, click Clear all on the selected items panel, or press Esc to clear the selection.

## Cell Groups

Each cell can belong to either one cell group or no cell groups. A cell cannot belong to multiple cell groups.

### To create a cell group:

- Select cells in the spatial view window, scatterplot, or heatmap. All selected cells are listed in the selected items panel. Cell indexes and identification numbers of the selected cells are also listed. The number of selected cells is listed in parentheses next to the panel title Selected items.
- Click on a cell in the selected items panel list and its location is identified by crosshairs in the spatial view window and scatterplot.
- To save selected cells as a cell group, enter a cell group name in the text box at the bottom of the selected items panel and click Save group. The cell group appears in the custom cell groups list in the spatial features panel. The number of cells in the group is listed in parentheses next to the cell group name.

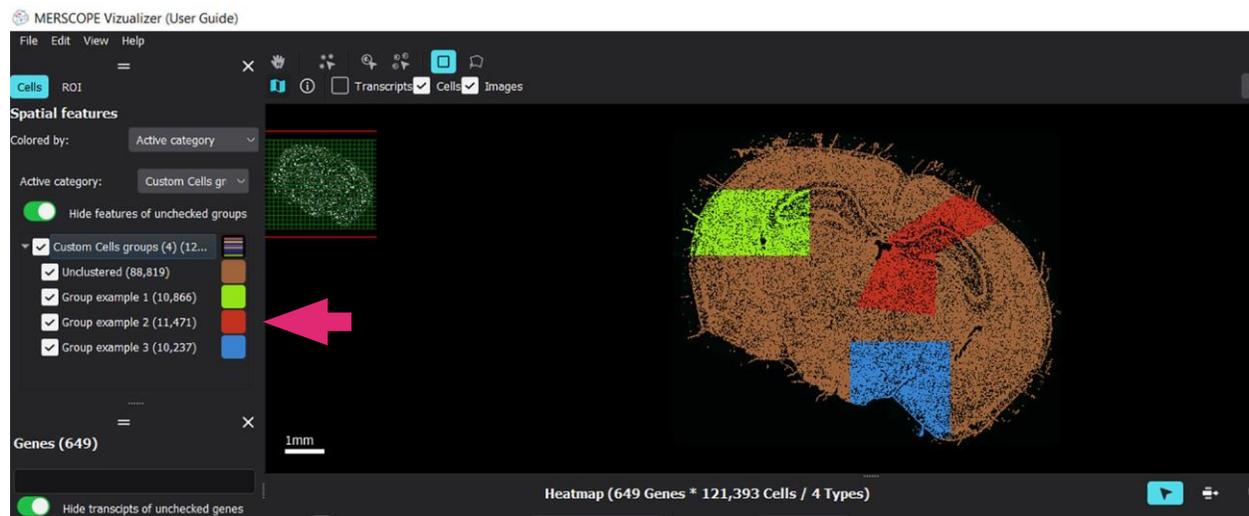

- *The newly created group always appears in the custom cell group category, regardless of which active category is selected.*
- *If a new cell group is created and the cell already belongs to a previously-defined cell group, the cell belongs to the newer group and is deleted from the previous group.*
- *After saving a cell group, the MERSCOPE Vizualizer software clears the selection.*
- *It is not possible to exclude (deselect) individual cells from a saved group.*

### To import a cell group:

- With an existing dataset open, export AnnData ([File > Export Data](#)) and export the cells feature set.

→ Import AnnData into a Jupyter notebook ([Open Custom Cell Clustering Analysis](#)).

```
In [1]: import scanpy as sc


In [13]: import pandas as pd
import numpy as np

In [4]: adata = sc.read_h5ad('../MsBrain_custom_cell_group.hdf5')
adata

Out[4]: AnnData object with n_obs × n_vars = 121393 × 649
obs: 'Custom Cells groups', 'Datasets', 'volume', 'center_x', 'center_y'
uns: 'Custom Cells groups_colors', 'Datasets_colors'
```

→ To create a new cell group, generate a new column of cell metadata.

– *In the example below, a new column of metadata (named new-group) was formed containing cells expressing the Slc17a7 gene at a high or low level.*

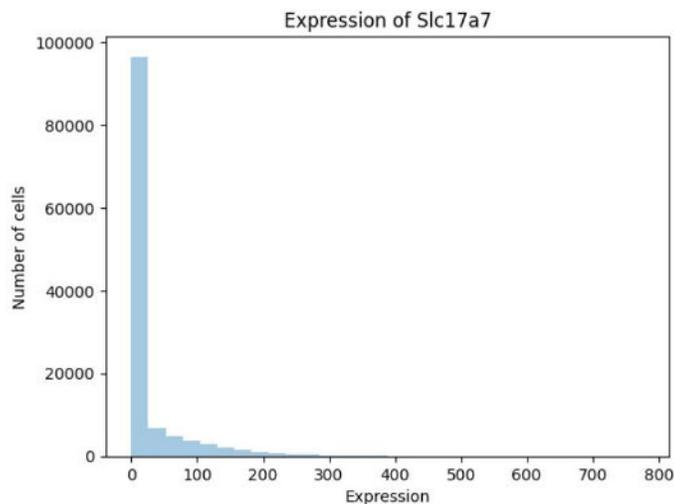

```
In [32]: ser_exp = adata.to_df()[gene_of_interest]
keep_cells = ser_exp[ser_exp > 30].index.tolist()

In [33]: adata.obs.loc[keep_cells, 'new-group'] = 'high'

In [34]: adata.obs['new-group'].value_counts()

Out[34]: new-group
low      98145
high     23248
Name: count, dtype: int64

In [35]: adata.write_h5ad('../MsBrain_new_group.hdf5')
```

→ The Jupyter notebook saves this AnnData to a new .HDF5 file.

→ Import this AnnData into the Vizualizer software ([File > Import Data](#)).

→ Select the new column of cell metadata under the list of categories (e.g., new-group in the example above).

→ The new cell category is available for selection and manipulation within the Vizualizer software.

### Checked Cell Groups

- Select an active category from the menu in the spatial view window. The corresponding cell group tree structure appears below in the window. Click the arrow to the left of the parent cell group to expand the tree structure.
- Checked cell groups appear in the spatial view window and scatterplot, and are also included in the heatmap.
- To select all cells in a checked cell group, click on the cell group name in the spatial features panel. Selected cells are shown in the spatial view window, scatterplot and heatmap, and are also listed in the details panel.
- In the case of multilevel hierarchy, checking a parent group applies the same status to all child groups.

### Unchecked Cell Groups

- Unchecked cells are hidden or colored gray in the spatial view window and scatterplot, and not included in the heatmap.
- Hide/unhide unchecked cell groups in the spatial view window and scatterplot by toggling the Hide features of unchecked groups slider in the spatial features panel.
- In the case of multilevel hierarchy, unchecking a parent group applies the same status to all child groups.

### Cell Group Settings

- To color any cell group, left-click the color spot to the right of the group name and select a color.
  - *The color is applied to the cell group in the spatial view window, scatterplot, and heatmap.*
- Right-click on any cell group name to:
  - Add/edit description of the cell group.
  - Reset color back to the default color.
  - Edit categories to rename or remove imported categories.
  - Remove group. The option to remove any custom cell group created in the MERSCOPE Vizualizer software. Default or imported cell groups do not have the remove option. Cells from the removed group return to the unclustered group.
  - Copy to clipboard. Copies the cell identification numbers in the cell group to clipboard. It is pasted as a comma delimited list.
  - Export. Exports the cell identification numbers to a .CSV file format.

→ Right-click on an active category name to:

- Reset color for all groups in the category back to the default colors.
- Edit categories to rename or remove imported categories.

### Cell Color Options

Cells can be colored according to cell category or a number of gene expression level characteristics.

Cells can be colored according to:

- Active category
- Count
- Z-score
- Normalized
- Normalized z-score

→ Cells are colored by active category by default. To change the coloring mode, select the option from the Color by menu at the top of the spatial features panel.

– *Active category applies the cell group colors defined by the color spot to the right of the cell group names.*

→ Selection of an expression level characteristic reveals a color scheme menu.

→ The three-color scheme based on minimum, center, and maximum values for each characteristic can be defined. Click on a color spot to change the color. Click on the center spot to convert to a two-color scheme (spot becomes transparent). Click again on the center spot to convert back to a three-color scheme and/or change the color.


→ Adjust the Brightness slider to define the minimum and maximum thresholds for the characteristic.

– *The expression level characteristics for every cell are calculated dynamically based on the total expression of the checked genes. To color based on the expression of a single gene or a subset/group of genes, uncheck all genes and then check a single gene or subset/group of genes.*

## SPATIAL FEATURES PANEL - REGIONS OF INTEREST

By default, there is one empty ROI feature set (named ROI) in the MERSCOPE Vizualizer software. ROI feature sets can be renamed or added via the [Edit > Manage ROI Sets](#) menu.

### Select an ROI Feature Set

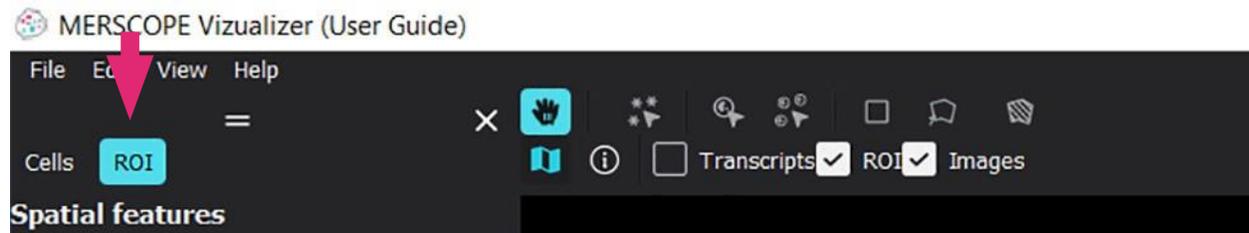


ROI feature sets and the cells feature set ([Select the Cells Feature Set](#)) are individually addressable by selection of a feature set at the top of the spatial features panel. The selected items panel also indicates the active feature set, i.e., the spatial feature set selected in the spatial features panel. The active feature set is always displayed in front of other feature sets in the spatial view window and selection tools are only applicable to the active feature set.

### Selection Toolbar (ROIs)

An ROI feature set layer in the spatial view window must be turned on to see ROIs. To toggle ROI feature set layers on and off, check/uncheck an ROI feature set layer in the [Spatial Settings](#) menu or in the shortlist of layers above the top left of the spatial view window (available when an ROI feature set is selected).

When an ROI feature set is selected in the spatial features panel, a Draw ROI polygon icon (6) appears in the selection toolbar above the top left of the spatial view window; the selection tools also work in the scatterplot.

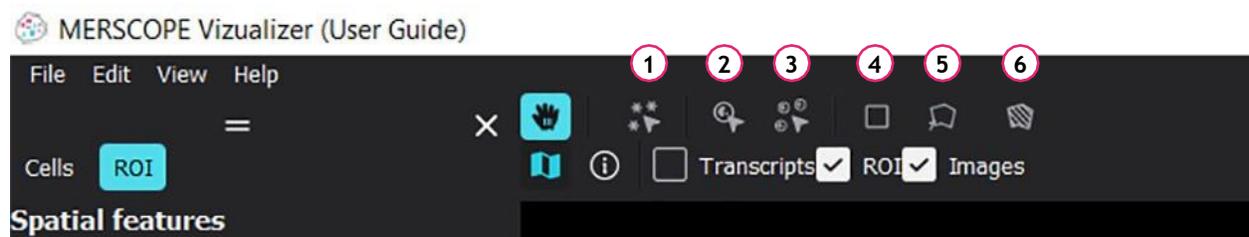

→ Multi-transcript selection (1) enables users to select all instances of a gene transcript at the same time. Click away from a selected transcript or press Esc and all [Checked Genes](#) transcripts reappear.

NOTE: Multi-transcript selection works on any LOD (Level of Details) for .VZG format, but for .VZG2 datasets multi-transcript selection works only on LOD 0.

→ Single-feature selection ([Name of ROI feature set]) (2) enables users to select a single ROI. Click/click again to select/deselect an individual ROI; click Clear all on the selected items panel, or press Esc to clear

the selection. Select multiple ROIs by Ctrl-clicking single ROIs one-by-one. To deselect a ROI from the selection, Ctrl-click in the spatial view window, scatterplot, or heatmap (as long as the Group by active category slider is toggled off in the [Heatmap Toolbar](#)).

- Multi-feature selection ([Name of ROI feature set]) (3) enables users to select all ROIs in an ROI group. Click/click again to select/deselect an ROI group; click Clear all on the selected items panel to clear the selection. Select multiple ROI groups by Ctrl-clicking ROI groups one-by-one.
- Rectangle selection (4). Drag and click to define a rectangle over the display to select ROIs at least partially within the rectangle. Press Ctrl and draw another rectangle to add ROIs to the previous selection. Draw another rectangle without Ctrl away from the original to reset the selection. Draw another rectangle within the first to select the subset of ROIs. Switch to the single-feature selection tool and click in a blank space in the spatial view window, click Clear all on the selected items panel, or press Esc to clear the selection.
- Polygonal lasso selection (5). Click each point to draw a polygon over the display. Close the polygon to select the ROIs at least partially within the polygon. Press Ctrl and draw another polygon to add ROIs to the previous selection. Draw another polygon without Ctrl away from the original to reset the selection. Draw another polygon within the first to select the subset of ROIs. Switch to the single-feature selection tool and click in a blank space in the spatial view window, click Clear all on the selected items panel, or press Esc to clear the selection.
- Draw ROI polygon (6). Click each point to draw a polygon over the display. Close the polygon and a dialog box appears.

### ROI Construction and ROI Groups

- To draw an ROI, use the Draw ROI Polygon tool. Click each point to draw a polygon over the display. Close the polygon and a dialog box appears. Press Esc before closing the polygon to reset the drawing.
- Enter a name for the newly drawn ROI.
- Select the group that the ROI should be added to or create a new group. To create a new group, overwrite the default group name in the ROI group box. To select from existing groups, click in the ROI group box and a list of existing ROI groups appears.
- Click Cancel to remove the ROI just drawn.
- *When an ROI feature set is selected in the spatial features panel, all ROIs in any selected group are listed in the [Selected Items Panel - ROI](#). The number of selected ROIs in the group is listed in parentheses next to panel title Selected items. Individual ROIs can be selected by clicking the ROI in the list and the center of the ROI is identified by crosshairs in the spatial view window.*

- *To save selected ROIs as a new ROI group, enter an ROI group name in the text box at the bottom of the selected items panel and click Save group. The ROI group appears in the custom ROI groups list in the spatial features panel. The number of ROIs in the group is listed in parentheses next to the ROI group name.*
- *To delete an ROI, right-click on the ROI name in the selected items panel.*
- *If a new ROI group is created and an ROI already belongs to a previously-defined ROI group, the ROI belongs to the newer group and is removed from the previous group. If the ROI group is empty (no ROIs) after moving an ROI, the group will disappear from the list in the spatial features panel.*
- *After saving an ROI group, the MERSCOPE Vizualizer software clears the selection.*
- *It is not possible to exclude (deselect) individual ROIs from a saved group.*

### Checked ROI Groups

- Checked ROIs appear in the spatial view window and scatterplot, and are included in the heatmap.
- To select all ROIs in a checked ROI group, click on the ROI group name in the spatial features panel. The ROIs in the group are listed in the selected items panel.
- In the case of multilevel hierarchy, checking a parent group applies the same status to all child groups.

### Unchecked ROI Groups

- Unchecked ROIs are hidden or colored gray in the spatial view window and scatterplot and not included in the heatmap.
- Hide/unhide unchecked ROI groups by toggling the Hide features of unchecked groups slider in the spatial features panel.
- In the case of multilevel hierarchy, unchecking a parent group applies the same status to all child groups.

### ROI Group Settings

- To color any ROI group, left-click the color spot to the right of the group name and select a color.
- *The color is applied to the ROI group in the spatial view window, scatterplot, and heatmap.*
- Right-click on any ROI group name to:
  - Add/edit description of the ROI group.

- Reset color back to the default color.
- Edit categories to rename or remove imported categories.
- Remove group. The option to remove any custom ROI group created in the MERSCOPE Vizualizer software. Default or imported ROI groups do not have the remove option. ROIs from the removed group return to the unclustered group.
- Copy to clipboard. Copies the ROI names in the ROI group to clipboard. It is pasted as a comma delimited list.
- Export. Exports the ROI names to a .CSV file format.

→ Right-click on an active category name to:

- Reset color for all groups in the category back to the default colors.
- Edit categories to rename or remove imported categories.

### ROI Color Options

ROIs can be colored according to ROI category or a number of gene expression level characteristics.

ROIs can be colored according to:

- Active category
- Count
- Z-score
- Normalized
- Normalized z-score

→ ROIs are colored by active category by default. To change the coloring mode, select the option from the Color by menu at the top of the spatial features panel.

– *Active category applies the ROI group colors defined by the color spot to the right of the ROI group names.*

→ Selection of an expression level characteristic reveals a color scheme menu.

→ The three-color scheme based on minimum, center, and maximum values for each characteristic can be defined. Click on a color spot to change the color. Click on the center spot to convert to a two-color scheme (spot becomes transparent). Click again on the center spot to convert back to a three-color scheme and/or change the color.

→ Adjust the Brightness slider to define the minimum and maximum thresholds for the characteristic.

– *The expression level characteristics for every ROI are calculated dynamically based on the total expression of the checked genes. To color based on the expression of a single gene or a subset/group of genes, uncheck all genes and then check a single gene or subset/group of genes.*

## SCATTERPLOT


The MERSCOPE Vizualizer software has a single scatterplot. The scatterplot can display two-dimensional arrays, such as cells in dimensionality reduced gene expression space (e.g., UMAP/ tSNE) or other user-imported two-dimensional arrays (e.g., cell volume versus transcript count). Clustering modes are imported with annotation data (AnnData). The active feature set is indicated in the scatterplot if features are colored by expression. If features are colored by active category, this category is shown in the scatterplot.. If there are no two-dimensional data for an ROI feature set, the scatterplot will be empty.

- *Only cells or ROIs from the experiment that have imported annotation are shown in the scatterplot. E.g., if some cells or ROIs were filtered during clustering, they are not displayed in the scatterplot.*

### Translate and Zoom

- In the zoom/pan cursor mode (hand icon) from above the spatial view window ([Navigation and Views](#)), click and drag the cursor in the scatterplot to translate the content.
- To zoom in/out of the content in the scatterplot, rotate the mouse wheel.

### Select Array

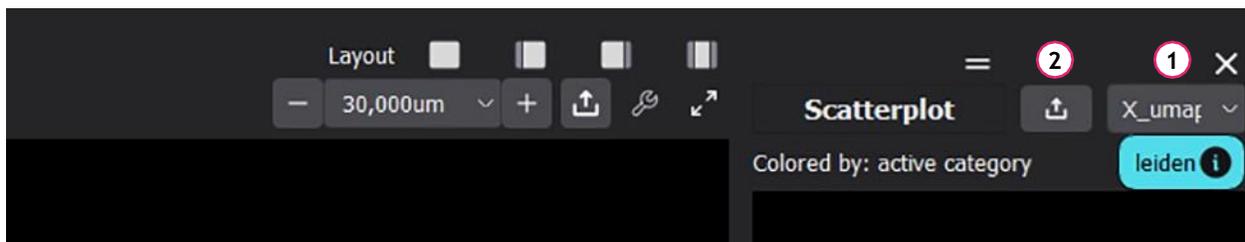

- If multiple arrays were imported, select the array visible in the panel from the menu above the scatterplot (1).

### Export Image

- Click the export icon (2) above the scatterplot to export a .png file of the content.
- Changing the resolution factor of the .png file changes the resolution and size of the image file. The range of resolution factors depends on the rendering capabilities of the computer and the actual size of the exported spatial view window (a smaller window can be exported with a higher resolution factor).

### Selection Toolbar (Cells or ROI)

Selection toolbar functions can be applied to the scatterplot in the same way as in the spatial view window. Refer to [Selection Toolbar \(Cells\)](#) or [Selection Toolbar \(ROIs\)](#) for more information.

Cell groups or ROI groups can be created in the scatterplot in the same way as in the spatial view window. Refer to [Cell Groups](#) or [ROI Construction and ROI Groups](#) for more information.

## GENES TREE

The genes panel controls the representation of gene products - transcripts - in the MERSCOPE Vizualizer software.

In the spatial view window, transcripts are represented by dots and all transcripts that are products of the same gene have the same color.

The transcripts layer in the spatial view window must be turned on to see transcripts. The transcripts layer is visible by default when an experiment is loaded. To toggle the transcripts layer on and off, check/uncheck the transcripts layer in the [Spatial Settings](#) menu or in the shortlist of layers above the top left of the spatial view window.

### Checked Genes

- All genes included in an experiment are listed in the genes panel. The number of genes in the gene panel is listed in parentheses next to the software panel title Genes and next to All in the list. The number of transcripts of each gene in the experiment (both inside and outside cell boundaries) is listed in parentheses next to the gene name. The number of genes in [Gene Groups](#) are listed in parentheses next to the gene group.
- To visualize all transcripts of a gene/gene group, check the box next to the gene/gene group name.
- Checked gene transcripts appear in the spatial view window.
- Checked genes are displayed in the heatmap.
- Selected genes are also highlighted in the heatmap ([Heatmap Highlighting](#)). Click on the gene name in the genes panel to select the gene.
- In the case of multilevel hierarchy, checking a parent group applies the same status to all child groups.

### Unchecked Genes

- Unchecked genes are hidden or colored gray in the spatial view window and scatterplot, and not included in the heatmap.
- Hide/unhide unchecked genes by toggling the Hide transcripts of unchecked genes slider in the genes panel.
- In the case of multilevel hierarchy, unchecking a parent group applies the same status to all child groups.

## Gene Color Settings

→ To change the color of gene transcripts, left-click the color spot to the right of the gene name and select a color.

## Gene Settings

→ Right-click on any gene name/color spot to see a menu to:

- Add/edit description of the gene.
- Move to another gene group - select the gene group from the menu or create a new group.
- Reset color back to the default color.
- Copy to clipboard. Copies the gene group, gene name, and color information in the gene to clipboard. It includes any description of the gene. It is pasted as a comma delimited list. The gene must be highlighted (left-click) prior to right-click to copy (otherwise information is not copied).
- Export. Exports the gene group, gene name, gene description, and color information to a .CSV file format. The gene must be highlighted (left-click) prior to right-click for export (otherwise the exported file will not contain information).
- Import. Gene groups, description, and color can be imported from a .CSV file. Click Import to upload a .CSV file, which must contain 4 columns: gene group, gene name, gene description, gene color; gene names are immutable. It is recommended to first export a .CSV file, edit it, and then import it.

## Gene Highlighting

- Left-click a gene or gene group to highlight it in the genes tree.
- Ctrl-click to highlight multiple genes.
- Right-click to move selected gene(s) to another group or create a new group.
- Highlighted genes are marked by a blue bar in the genes tree.
- Use the up/down arrow keys to move highlighting up/down in the genes tree.
- Unchecked genes are not highlighted if Hide transcripts of unchecked genes is toggled on.
- In the spatial view window:
  - Highlighted gene(s) are colored.
  - Non-highlighted genes are gray.
  - Hide/unhide selected genes by toggling the Hide Transcripts of Unselected Genes slider in the Genes panel.
- In the heatmap:
  - Highlighted gene(s) are marked by a blue bar ([Heatmap Highlighting](#)).
- A gene must be checked in order to be shown in the visualization regardless of highlighting.


## Gene Groups

Each gene can belong to either one gene group or no gene groups. A gene cannot belong to multiple gene groups.

Genes without a group are designated ungrouped. A gene group can be created from checked genes, from highlighted genes, or an empty gene group can be created. If a gene group is deleted, the genes in the group automatically return to the ungrouped status. It is not possible to delete a gene from the genes panel.


To create a gene group from checked genes:

- Check genes in the genes panel.
- To save checked genes as a gene group, enter a gene group name in the text box at the bottom of the genes panel and click Save. The gene group appears in the tree structure in the genes panel. The number of genes in the group is listed in parentheses next to the gene group name. The number of transcripts of each gene in the experiment (both inside and outside cell boundaries) is listed in parentheses next to the gene name.

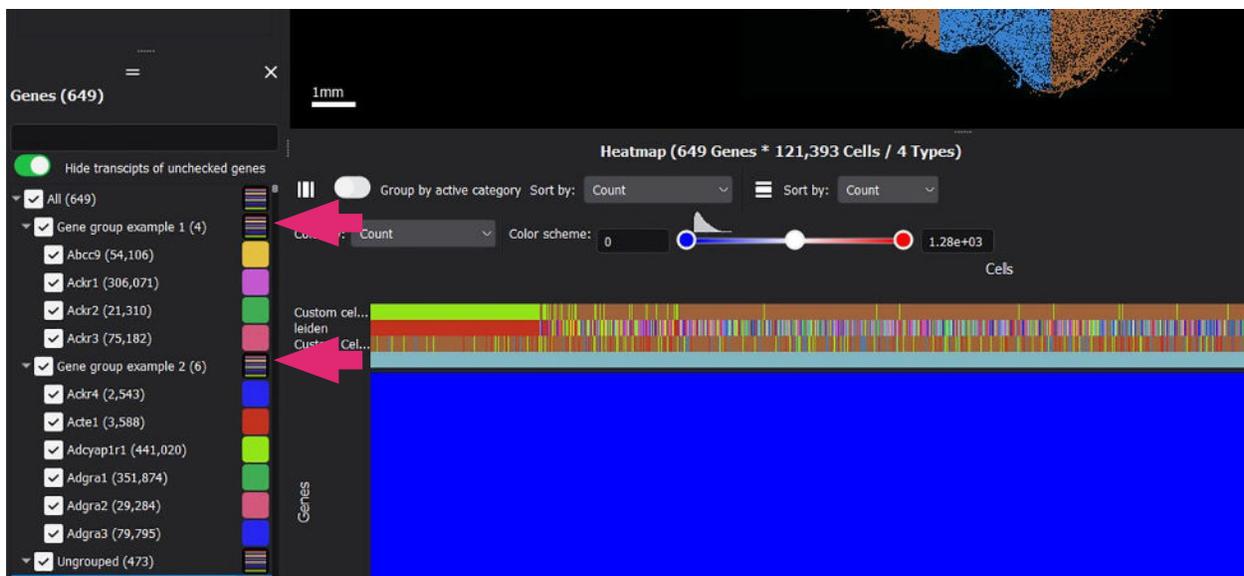

- Click the arrow to the left of the parent gene group to expand the tree structure.
- *If a new gene group is created and the gene already belongs to a previously-defined gene group, the gene belongs to the newer group and is deleted from the previous group.*
- To create an empty gene group, ensure no genes are selected, enter a gene group name in the text box at the bottom of the genes panel and click Save.
- *An empty gene group can be useful if users want to create a gene group in a stepwise fashion, adding one gene at a time, for example.*
- Gene groups can be created by right-click from highlighted genes (Move to > Create new

group).

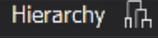- Gene groups can be created from the heatmap by clicking on an individual cluster or by clicking the hierarchy button  in the genes panel when hierarchical clustering is selected in heatmap toolbar.
- Gene groups can be highlighted in the heatmap.

### Search for Genes

- Use the search bar at the top of the genes panel to search for genes in all groups.


### Gene Group Settings

- Right-click on any gene group name/color spot to see a gene group menu to:
  - Color by group(s). Changes all genes in the same group to the same color. To change the color of a gene group, left-click the color spot to the right of the gene group name and select a color.
  - Reset color back to the default color.
  - Delete group. Provides the option to delete any gene group created in the MERSCOPE Vizualizer software. When a gene group is deleted, the genes in the group return to the ungrouped status.
  - Copy to clipboard. Copies the gene group, gene name, and color information of all genes in the gene group to clipboard. It is pasted as a comma delimited list. The gene group must be highlighted (left-click) prior to right-click to copy (otherwise information is not copied).
  - Export. Exports the gene group, name, gene description, and color information to a .CSV file format. It is possible to save the different types of genes grouping, e.g., grouping based on hierarchical clustering, DEG analyses, etc. The gene group must be highlighted (left-click) prior to right-click for export (otherwise the exported file will not contain information).
  - Import. Gene groups, description, and color can be imported from a .CSV file. Click Import to upload a .CSV file, which must contain 4 columns: gene group, gene name, gene description, gene color; gene names are immutable. All genes from gene panel will be regrouped according to the information in the .CSV file. It is recommended to first export a .CSV file, edit it, and then import it.

## HEATMAP


The MERSCOPE Vizualizer software has a single heatmap, where users can view gene expression data according to spatial feature categories and groups. All cells from all datasets are shown in the heatmap and scatterplot. To view cells from a specific dataset(s) in an experiment assembled from multiple datasets, select Datasets in the [Active Category](#) menu in the spatial features panel and uncheck any listed dataset to hide the cells from that dataset.

### Navigation and Selection

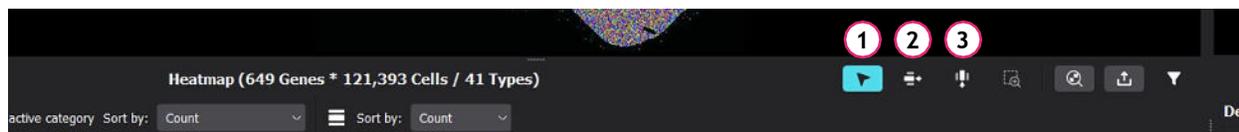


- In the zoom/pan cursor mode (arrow icon) (1), click and drag the hand cursor in the heatmap to translate the content of the window.
- In the gene selection cursor mode (row icon) (2), click a gene in the heatmap to select it. Information on the selected gene appears in the details panel.
- In the feature selection cursor mode (column icon) (3), click a feature in the heatmap to select it. Information on the selected feature appears in the selected items panel.

### Zoom and Scale

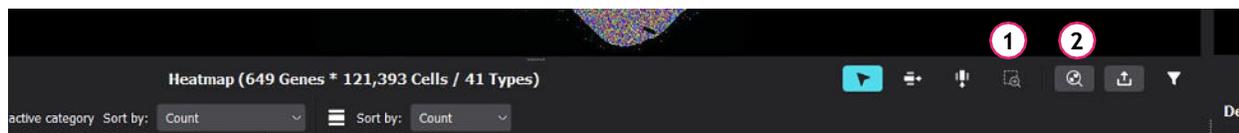

- To zoom in/out of the content in the heatmap, rotate the mouse wheel.
- In the zoom to rectangle mode (1), draw a rectangle over the heatmap to expand the contents of the rectangle.
- Use the reset zoom mode (2) to visualize the entire heatmap.


### Export

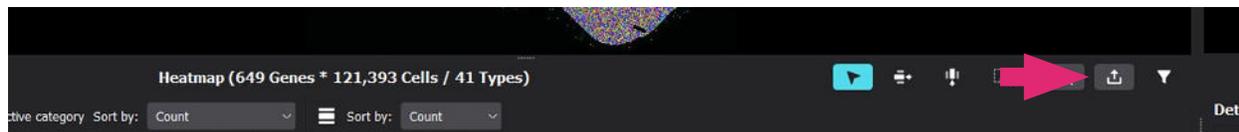

- Click the export icon to export a .png file of the heatmap or export the filtered expression matrix to a .CSV file or .PARQUET file.

## Heatmap Toolbar

The spatial feature set (cells or ROI) selected in the spatial features panel is visible in the heatmap and the data are ordinarily displayed per feature. Only [Checked Cell Groups](#) or [Checked ROI Groups](#), and [Checked Genes](#) are visible in the heatmap.

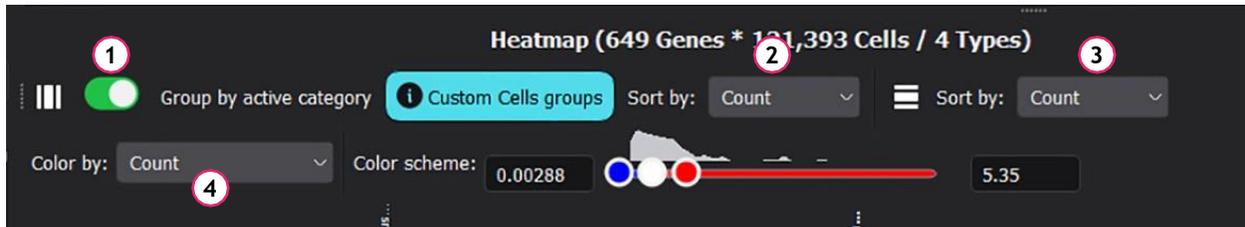

- The active category selected in the cells/ROI feature set is displayed in the heatmap. To display data according to cell/ROI groups in a category, toggle on the Group by active category slider (1).
- All cell categories are shown in the heatmap bar (at the top) in the order they appear in the active category list in the spatial features panel. When the Group by active category slider (1) is toggled on, only the active category is shown in the heatmap bar.
- Cells/ROIs can be sorted according to count or by count within each category (count+category) via the feature sort by menu (when the cells feature set is selected) (2).
- Genes can be sorted according to count, variance, or hierarchical clustering via the gene sort by menu (3). If hierarchical clustering is selected, it is possible to create gene groups via the hierarchy button in the genes panel ([Gene Groups](#)).
- When the Group by active category slider (1) is toggled on, the feature sort by menu (2) enables sorting by hierarchical clustering or count. The sliders on each axis of the heatmap (bottom left for genes and top right for category) adjust the level of the hierarchical dendrograms that are displayed as brackets on either axis (sliders are only visible when hierarchical sorting is selected).
- If selected genes or feature groups are changed, orange warning triangles  appear next to the sort by menus and the dendrograms appear orange. Click the recalculate hierarchical clustering button to the top left of the heatmap (indicated by an orange warning triangle/circular arrow) to adjust the clustering according to the new selections.
- Spatial features (cells or ROI) in the heatmap can be colored according to gene expression level. Spatial features can be colored according to count, Z-score, normalized, and normalized z-score (4).
- The three-color scheme based on minimum, center, and maximum values for each characteristic can be defined. Click on a color spot to change the color. Click on the center spot to convert to a two-color scheme (spot becomes transparent). Click again on the center spot to convert back to a three-color scheme and/or change the color.

→ Adjust the slider to define the minimum and maximum thresholds for the characteristic.

## Heatmap Settings

To manage properties of the heatmap, toggle the heatmap settings menu on and off via the funnel icon above the top right of the heatmap.



→ Adjust the Variance of highly variable genes slider to set a cutoff for genes to be displayed in the heatmap, and click Apply. When adjusting the slider, the number of kept genes appears under the box, next to an information icon.

– *After clicking Apply, the genes below cutoff are unchecked in the genes panel. To reset the genes included in the heatmap, re-check the genes in the genes panel.*


→ Adjust the Z-score of differentially expressed genes (DEG) slider to set a cutoff for genes to be displayed in the heatmap, and click Apply. When adjusting the slider, the number of kept genes appears under the box next to an information icon.

- Only checked gene groups are taken into account in DEG calculation.
- At least two cell groups or ROI groups should be checked for Z-score or DEG calculation.
- DEG calculation for a large number of groups may take some time (up to 60 sec).

– *If the Z-score of differentially expressed genes slider is not visible, click Recalculate DEG and the slider will appear. This may take some time.*

– *After clicking Apply, the genes below cutoff are unchecked in the genes panel. To reset the genes included in the heatmap, re-check the genes in the genes panel.*

## Heatmap Highlighting

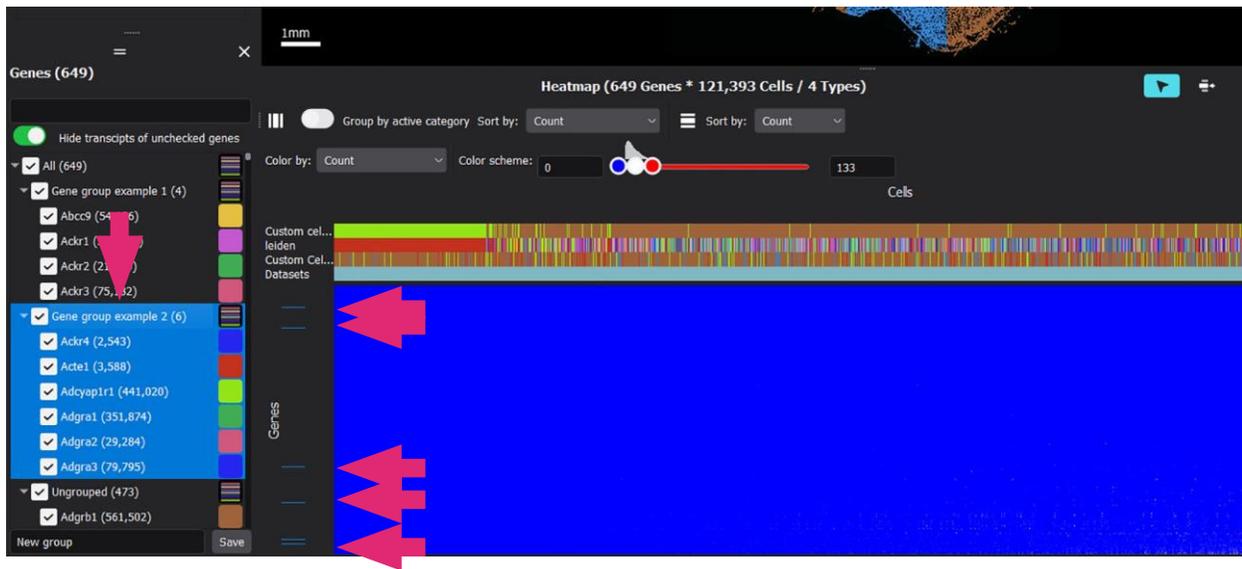

Locations of selected and checked items are highlighted in the heatmap by a blue bar on the applicable axis. Items must be checked to be highlighted (gene group member highlighting illustrated). Also refer to [Gene Highlighting](#).

Highlighting on the y-axis (genes):

- Individual genes are highlighted when a gene transcript is selected in the spatial view window.
- Individual genes are highlighted when a gene is selected in the genes panel.
- Gene group members are highlighted when a gene group is selected in the genes panel.
- Individual genes are highlighted when a gene name (label) on the left side of the heatmap is selected.
- Use Ctrl-click to select multiple genes.

Highlighting on the x-axis (cells or ROIs):

- Individual cells or ROIs are highlighted when a cell or ROI name (label) at the top of the heatmap is selected. Use Ctrl-click to select multiple individual cells or ROIs. The selected (and highlighted) cells or ROIs are listed in the selected items panel.
- Individual cells or ROIs are also highlighted when a cell or ROI is selected in the spatial view window or scatterplot.
- Cell or ROI group members are highlighted when selected in the spatial features panel, or in the spatial view window or scatterplot using the multi-feature selection tool.

Cell/ROI names and custom cell names can be copied from heatmap axes by right-click on the name.